

TWO POSITION MICRO-ELECTRIC VALVE ACTUATOR

Technical Note 421

Control Module Serial Numbers Starting with E2CA*

The two position micro-electric actuator consists of a control module, a stepper motor/gearbox assembly, a manual controller (use is optional), a universal AC input (100-240 VAC, 50-60 Hz) to 24 VDC power supply, and the interconnecting cables. The actuator is for use with all Valco and Cheminert two position valves with stops, toggling them back and forth between Positions A and B and outputting a position indication at the end of each move. (Position A is the counterclockwise position as seen when facing the output shaft of the motor; Position B is clockwise.)

The actuator is self-adjusting—no manual valve alignment is necessary. The end points of the valve travel are automatically located during a user-prompted initialization sequence which must be run each time a valve is mounted on the motor assembly (refer to page 4).

Actuator torque and speed are functions of the particular gearbox/motor/controller combination, which is designated as Model EQ, EH, EP, ED, ET, or EB in order of increasing torque/ decreasing speed. Models EQ, EH, ED, and ET have unique gearboxes; Models EP and EB have EH and ET gearboxes, respectively, with the motor traveling at a slower speed to provide increased torque. Model gear ratios vary from 3.14 to 22.56, capable of continuous operation of valves with turning torque requirements from 2.7 to 50 lb/in (0.3 to 5.7 Nm), as indicated in **Table 1**.

VALVE SWITCHING TIMES

The actual time required to switch the valve from Position A to Position B (Table 2) depends on two factors: the model number (reflecting the actuator speed) and the degree of rotation between positions.

*Serial numbers are on the underside of the control module. (Figure 2) Actuators of this series are further identified by a plastic control module housing. If your actuator control module has a metal housing and a serial number starting with EM2C, please refer to Technical Note 413.


Table 1: Recommended maximum loads

ACTUATOR MODEL	MAX RECOMMENDED VALVE TORQUE LOAD
EQ	2.7 lb/in (0.3 Nm)
EH	7.0 lb/in (0.8 Nm)
EP	9.0 lb/in (1.0 Nm)
ED	22 lb/in (2.5 Nm)
ET	40 lb/in (4.5 Nm)
EB	50 lb/in (5.7 Nm)

Table 2: Valve switching times in milliseconds

MODEL	NUMBER OF PORTS IN VALVE					
	4	6	8	10	12	14
EQ	105	75	70	65	55	50
EH	145	105	85	70	65	70
EP	220	125	110	90	75	65
ED	425	290	230	200	170	155
ET	650	450	360	300	265	240
EB	1500	1050	830	700	615	570

Figure 1: Actuator and controller connections

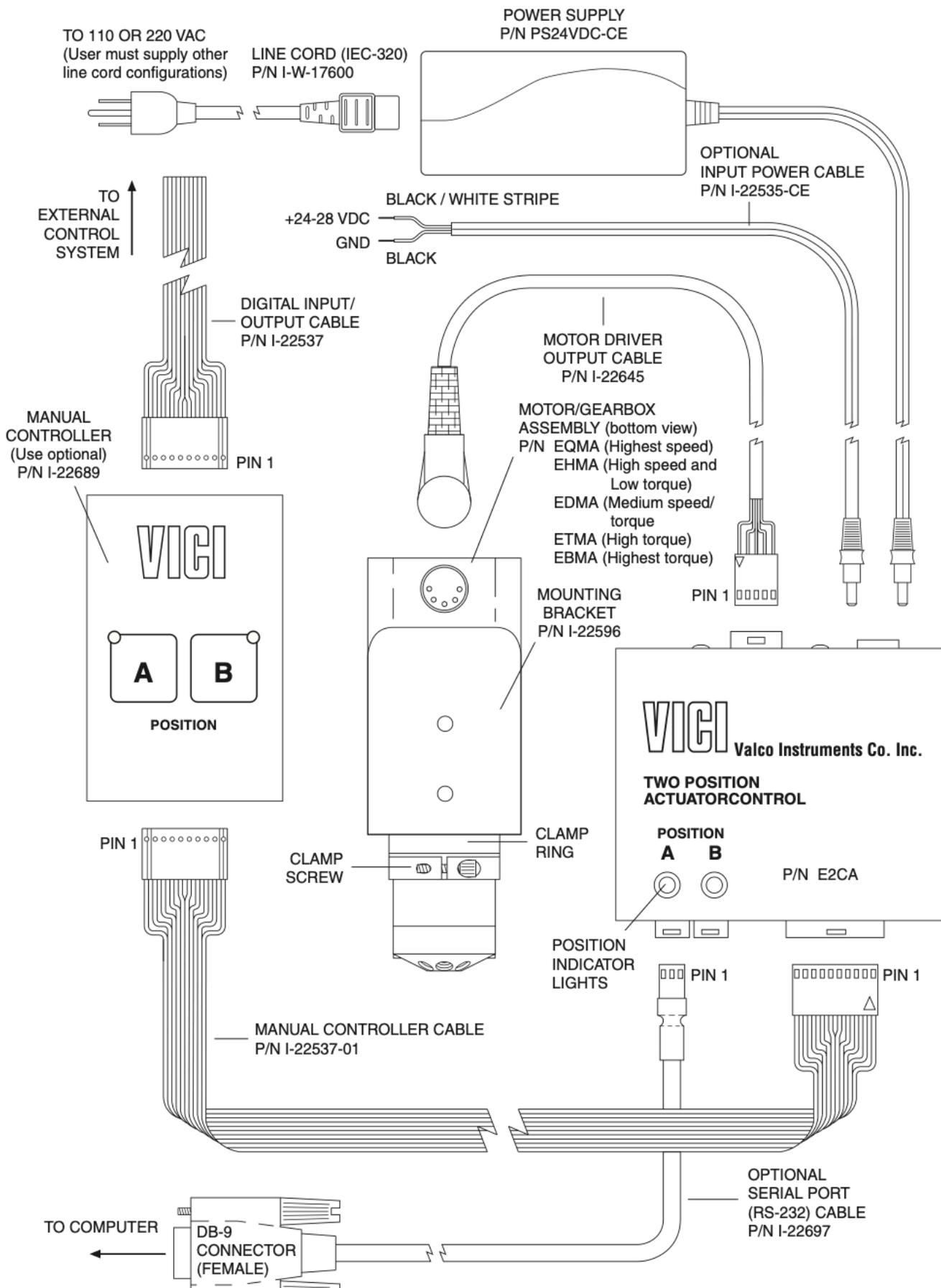

INSTALLATION AND USE

Getting Started

Figure 1 shows how to connect the various actuator components. There are four connectors on the control module, keyed and sized to prevent incorrect connection.

Cable and Connector Functions

Input power (24-28 VDC regulated) is supplied through a 2.5mm DC power jack; the inner pin is + voltage and the outer pin is ground. The average DC current requirement is 2.5 amps. Standby current draw is 60 milliamps. The actuator should not share a power supply with other noise-sensitive devices, as the high current draw may cause problems for these devices.

Motor driver output is through the five-pin connector: pins 1, 2, 4, and 5 carry the stepper motor phase drive signals. Pin 3 is grounded and tied to the cable shield to reduce electrical noise.

Two three-pin connectors are used for the serial port control. They are connected in parallel to facilitate implementation of multidrop applications. (Information on serial control of the actuator begins on page 4.)

The ten-pin connector is for digital inputs and outputs. A digital control system can operate and monitor the actuator through this port. The manual controller can be attached to the same connector and has a pass-through connection so it can be used along with an external control system.

Table 3

Pin #	Signal Description
INPUT POWER CABLE	
inner	+24 -28 VDC power
outer	Ground
MOTOR DRIVER OUTPUT CABLE	
1	Phase B
2	Phase B
3	Ground
4	Phase A
5	Phase A

Mounting

The actuator should be oriented so that any potential leakage of liquid from the valve or fittings flows away from rather than into the actuator (right). Figure 2 provides the mounting dimensions for the stepper motor/gearbox assembly and control module.


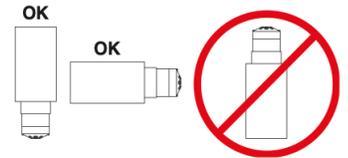


Figure 2: Stepper motor/gearbox assembly and control module

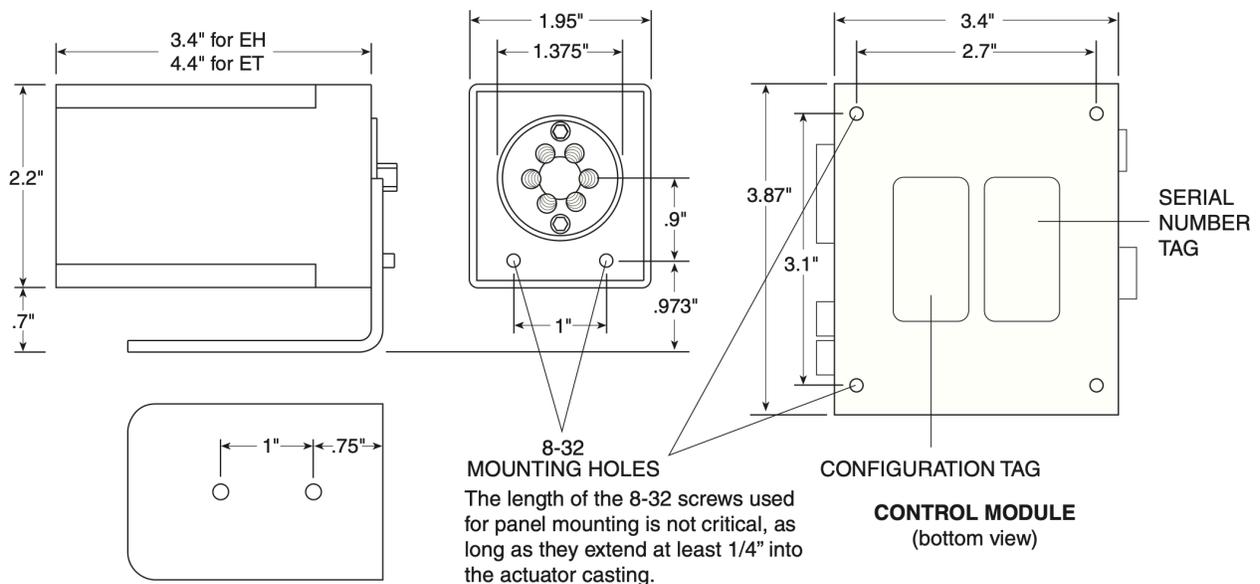

Initialization

Initialization is the procedure by which the actuator “learns” the stroke of the installed valve and its hardware. Any time a valve is removed and reinstalled, the actuator must be initialized by following these steps:

Hardware method:

1. Cycle the actuator twice with no valve, or with either end of the motor driver output cable unplugged.
2. Plug the cable back in, or put the valve back in the clamp ring. Orient the valve as desired and tighten the clamp screw.

Software method:

1. Send **IN**<enter> via serial communication. (Discussion of serial control begins below.)


Once you have started an initialization sequence by either method, the actuator will move at half its normal speed while it seeks the valve stops. Once it determines the proper travel, the actuator will return to normal speed. Cycle the actuator back and forth for four or five cycles, or until you hear the change in speed. If you don't hear the change, make sure that the clamp screw is tight.

Since the actuator checks the end points on each move, it will detect any inconsistencies. Positioning inconsistencies or errors are indicated when position outputs are deasserted or turned off (discussion of digital control begins on page 8).

SERIAL CONTROL OF THE ACTUATOR

Establishing Serial Communications

NOTE: For RS-485 information, refer to “Serial Control: RS-485 Option” on page 6.

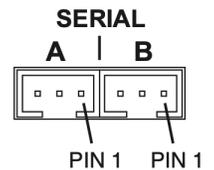

Items required:

- Terminal emulation or communication software such as Valco's VCom or Windows® Hyper- Terminal, running on a PC-compatible computer
 - Valco cable assembly I-22697 or equivalent
1. Connect the I-22697 cable to the actuator as indicated in **Figure 1**, and set the serial port at 9600 baud, no parity, 8 data bits, 1 stop bit, no hardware or software handshaking. Note: Valco's VCom software automatically sets the serial port to these default settings.)

Table 4

PIN #	RS-232	RS-485	DB9*
1	Ground	Ground	5
2	Transmit to host	A (-)	2
3	Receive from host	B (+)	3

2. With the software running, check the bi-directional communication link between the keyboard/ monitor of the computer and the serial port by typing **VR**<enter>. If the link is functioning and an actuator ID has not been set, a message similar to the following will appear on your monitor, giving the program number and date of the actuator firmware.

I-PD-E(t)F56R(xx) (t = actuator Model Q, H, P, D, T, or B; xx = revision letter and suffix)

08/06/2009 (Month first; zeros are included in the response)

If there is no response, it is possible that the ID has already been set. To force a response from a device with an unknown ID, type ***VR**<enter>. The asterisk is a wild card ID substitute which will elicit a response from all devices on line, no matter what their ID is.

Protocol

The serial protocol is an ASCII character string protocol. All commands are terminated with the carriage return (OD hex or '\r') character, which defines the end of commands for both inputs and outputs. Line feed characters (OA hex) sent to the device are ignored. When using a terminal program for communications, pressing the <enter> key will generate the carriage return character. Software flow control (Xon/Xoff) and hardware handshaking are not supported.

Table 5

SERIAL COMMANDS	
VR <enter>	Displays the part number and date of the firmware
GO(l) <enter>	Sends the actuator to Position <i>l</i> , where <i>l</i> is A or B
CW <enter>	Sends the actuator to Position A
CC <enter>	Sends the actuator to Position B
TO <enter>	Toggles the actuator to the opposite position
TT <enter>	Toggles the actuator to the opposite position, waits a preset delay time, then rotates back to the original position.
CP <enter>	Displays the current actuator position
DT(n) <enter>	Sets or displays the delay time, where <i>n</i> =0 to 65,535 milliseconds
ID(n) <enter>	Sets or displays the device ID to value <i>n</i> , from 0 to 9
ID(l) <enter>	Sets or displays the device ID to value <i>l</i> , from A to Z (RS-485 only)
SB(n) <enter>	Sets or displays the baud rate, where <i>n</i> = 2400 to 115,220 (default is 9600)
SM(n) <enter>	Sets or displays the digital input mode, where <i>n</i> = 1 to 4
SO(n) <enter>	Sets or displays the delay time before the position outputs are turned off, set in milliseconds to the closest 5 ms interval, from 0 to 30,000 ms. The outputs are always on (SO=0) by default.
CNT(n) <enter>	Sets or displays the move count, from 0 to 65,000
GV	Displays the measured DC input voltage
GVD	Displays the measured DC input voltage drop while moving
IN <enter>	Starts a re-initialization sequence
/? <enter>	Displays list of valid commands

Table 5 (previous page) includes commands to reposition the valve, to configure input and output functions, to set baud rates and other communication options, and to monitor the power status and the number of actuation cycles. All values that are set by serial port commands, such as **DT**, **ID**, **SM**, etc., are retained through power on/off cycles. Refer to special instructions for permanently setting the baud rate on page 6.

NOTE: Characters shown in parentheses indicate a user-defined number (*n*) or letter (*l*). If the command is issued without that character, the response displays the current setting; supplying the number or letter sets the actuator accordingly.

Using the Device ID Feature

Actuators are shipped from the factory in the RS-232 mode, with the ID feature disabled. When an ID is set, the actuator responds only to commands which begin with the correct ID prefix, and its transmit output is disabled when not in use. This allows up to 10 actuators to be controlled from one computer RS-232 port. A single command can be broadcast to all actuators by using an asterisk (*) as the command prefix. NOTE: Any broadcast command which elicits a response from the serial port (such as *VR or *ID) will elicit a combined and unintelligible response.

For permanent multidrop applications, the RS-485 mode (below) is the preferred solution. However, just as RS-232 control requires the host to have an RS-232 serial port, the PC host or control system must have an RS-485 port to communicate with the actuator in the RS-485 mode. Plug-in PCI cards with RS-485 ports or adapters that change an RS-232 signal to an RS-485 signal are available from several common electronic manufacturers. If your computer lacks a serial port, adapters which convert USB ports to RS-232 or to RS-485 are also readily available.

To set the ID of an actuator, connect it to a serial port as shown in Figure 1 on page 2.

Caution: When installing or replacing actuators on a shared serial port, make sure that no two devices have been set to the same ID number.


1. Remove all of the actuators from the serial daisy chain except the one for which you are setting the ID.
2. Type VR <enter>. You should get a response giving the firmware version, indicating that serial communication with the actuator is established. If there is no response, type *VR<enter> to see if the ID is already set. If there is still no response, check the cabling and connections.
3. To set an ID, type IDn<enter>, where *n* is the new ID, from 0 to 9.
To change an ID, type iIDn<enter>, where *i* is the current ID and *n* is the new ID.
To disable the ID feature, type iID*<enter>, where *i* is the current ID.

Setting a New Baud Rate

To permanently set a new baud rate for the actuator:

1. Establish communications with the actuator at the current baud rate.
2. Issue the command SBn to temporarily change the baud rate to the desired rate. If the power goes down at this point, the baud rate will revert to the last permanent setting.
3. Change the host computer to the same baud rate just set in the actuator, and verify that you can establish communication.
4. Re-issue the same SBn command as in Step 2, and the current baud rate will be made permanent.

Figure 3: Multiple actuators controlled by one computer via RS-232 or RS-485 (preferred)

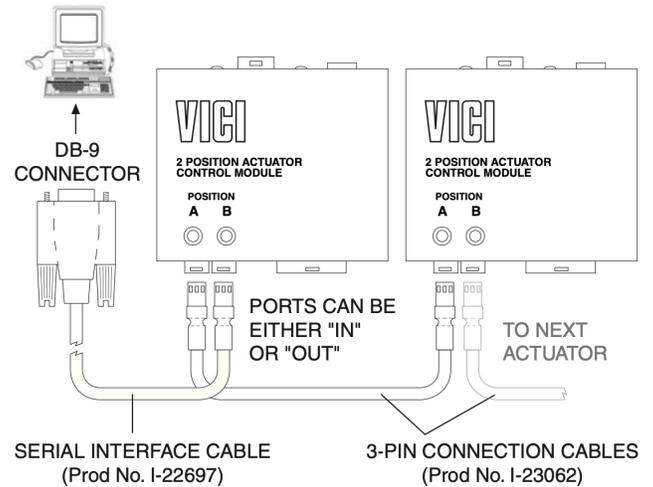

SERIAL CONTROL: RS-485 OPTION

When the serial port configuration switch is set to the "485" position, the serial port output is changed to the RS-485 mode. This also changes the software protocol, as described in the section entitled "Software" on the next page. You may wish to modify the configuration tag on the bottom of the control module to reflect this Comm change.


Setting the Serial Port Configuration Switch

1. Grip the lower section of the control module enclosure on the two opposing sides that have the curved release tabs.
2. Carefully depress the tabs while lifting the top section with the other hand. We recommend that the enclosure be opened in a static free environment following all proper ESD protection techniques.

NOTE: Once the control module cover has been removed, the printed circuit board can fall out.

3. Locate the slide switch marked "232 <- -> 485", and set it to 485. The slide switch labeled "Termination On <- -> Off" is usually left in the Off position, unless the wiring from the host control to the device is very long, and it is the last device at the end of a signal chain of devices.

Figure 4: Switch Locations

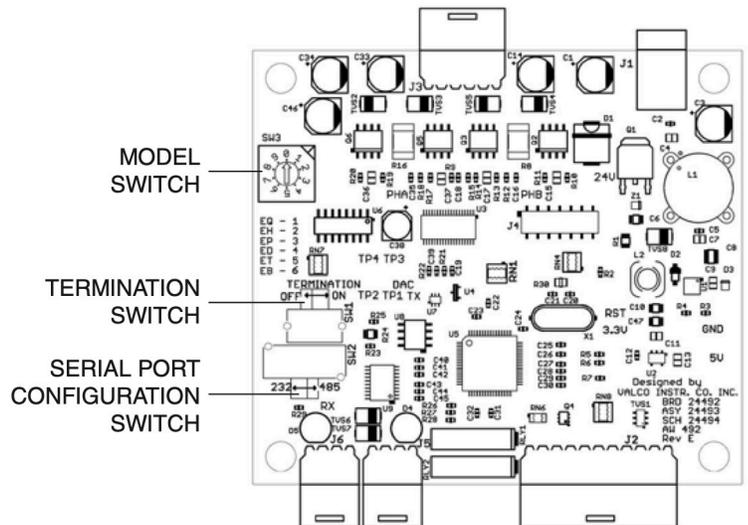

Caution: The rotary switch marked SW3 reflects which motor assembly is used, and should usually not be changed unless the control module is being dedicated to a different motor/gearbox assembly. If that is the case, refer to the section entitled "Setting the Model Switch" on page 9.

Hardware

The Valco implementation of RS-485 is half-duplex two wire with ground, and uses the same 3 pin output connector as RS-232. Pin 1 is ground, and pins 2 and 3 are signals 'A' (-) and 'B' (+), respectively.

RS-485 Cable

The cable should be 24 AWG twisted-pair with low shunt capacitance.

Configuring RS-485 Adapters

While each adapter is different, there are three issues to consider. The first consideration is two wire versus four wire application. Some adapters may have the capacity for both types of operation, and you may have to do some special wiring to support the two wire mode required for the Valco E2CA.

Another issue is related to flow control. Some adapters require the use of the serial port signal RTS to switch the adapter's RS-485 output from receive to transmit. In that case, you will need to enable the hardware handshaking control for that adapter to work.

The third issue to consider is termination. While most applications will work fine without any termination, if the cable length is hundreds of feet and the baud rate is set very high, then the termination is recommend for the

two farthest end points of the cable drop. It should not be used for devices in between. The actuator includes a switch which will enable and disable termination. (Refer to "Setting the Serial Port Configuration Switch" on the previous page.)

For details covering the implementation of your specific RS-485 hardware port or adapter, please consult the device's user manual.

Software

The RS-485 option involves three minor software adaptations to the RS-232 protocol. First, the ID range is extended to include the characters A through Z. (Upper and lower cases are treated as the same ID.)

The second change is that an ID is required (either numbers from 0 to 9 or letters from A to Z), and therefore must be included in all commands. If the actuator has already been set to an ID other than the factory default of **Z**, that ID will be retained; otherwise, the ID will be set to the **Z** default.

The third adaptation is that all commands must include a forward slash [/] as the start-of-message character. Example: The **VR**<enter> command used in the RS-232 protocol becomes **/ZVR**<enter> in the RS-485 protocol, where **Z** is the current ID of the device.

The RS-485 port on the host computer or controlling device generally includes terminating resistors, so only the actuator on the end of a long communication network needs to have its termination switch set to the On position.

DIGITAL CONTROL OF THE ACTUATOR

The digital input/output interface operates at 5 VDC signal levels and provides a limited amount of power that can be used for an external user interface. The asserted, or "ON", state for all output or input signals is a low voltage logic level, typically between 0 and .8 VDC. The deasserted, or "OFF", state is a high voltage logic level, typically between 4 and 5 VDC. This type of logic is referred to as "negative true" logic. All inputs have 5K ohm pull up resistors for a disconnected default "OFF" state.

Power Output

Pins 1 and 2 provide ground and +5 VDC, respectively, with a maximum allowable current draw of 100ma.

Caution: Do not connect the +5 VDC output to the power output on any external device, as this may cause permanent damage to the controller. Also be sure that anything connected to the +5 VDC connection does not draw more than 100 ma of current. Excessive current draw can cause the controller to malfunction.

Digital Outputs

There are two separate outputs for each position indication – a digital output and an isolated relay contact output for both Position A and Position B. Only one set of these outputs is asserted at a time. For a "Position A" indication, pin 3 is driven to a logic low and pins 7 and 8 are shorted for the isolated relay contact output, while pin 4 is at a logic high and pins 9

DIGITAL I/O CABLE	
Pin #	Signal Description
1	Ground
2	+5 VDC
3	Position A output
4	Position B output
5	Position A input
6	Position B input
7	Position A relay contact output
8	Position A relay contact output
9	Position B relay contact output
10	Position B relay contact output

and 10 are open. For a "Position B" indication, pin 4 is driven to a logic low and pins 9 and 10 are shorted while pin 3 is at a logic high level and pins 7 and 8 are open. If there is a positioning error due to valve sticking, clamp ring slippage, or re-initialization, both digital outputs are set to the "OFF" state (high logic level), and both relay contacts are open.

Digital Inputs

There are two digital inputs, pins 5 and 6, which can be driven by a 5 volt 1.5 milliamps TTL/CMOS logic signal, by an open collector driver, or by a switch or relay used to short the signal to the ground on Pin 1. All input signals must be asserted or deasserted for a minimum of 30 milliseconds to be properly recognized and processed. Any signal once asserted that is held or reasserted is ignored.

Caution: Do not connect any voltage signal above +5 VDC to these inputs as permanent damage to the device may result.

Digital Input Modes

Four input mode options are provided to expand the control flexibility of the actuator. The factory default setting is Mode 1. All four modes are described below. The mode can be set via the serial port using the **SMn** command and is retained even when the power is cycled to the controller. See the serial commands section on page 4-5 for details. NOTE: Changing the input mode does not affect the digital output functions.

Mode 1 (dual signal control mode)

In mode 1 (default) the digital inputs use the two input signals for position control. Asserting input pin 5 causes the actuator to go to Position A, and asserting input pin 6 sends it to Position B. If both signals are asserted at the same time, they will be ignored.

Mode 2 (single signal toggle mode)

Operation in this mode uses a single signal for position control with two options. Asserting pin 5 causes the actuator to toggle from the current to the opposite position. Asserting pin 6 causes the actuator to toggle to the opposite position, delay for a user settable period of time (the default is 100 ms), then toggle back to the original position. The delay time is set using the **DTn** serial port command. See the serial commands section on page 4-5 for details.

Mode 3 (state mode with enable)

This mode uses the pin 5 input as a state input, meaning that a low signal (asserted) on pin 5 causes the actuator to move to Position A. A high signal (deasserted) on pin 5 causes the actuator to move to Position B. Pin 6 is used as an enable/disable pin to prevent startup irregularities. The actuator is normally disabled when pin 6 is in its high (deasserted) state; it must be pulled low (asserted) to enable the actuator to move.

Mode 4 (state mode with enable)

This mode differs from Mode 3 only in the status of pin 6. As in Mode 3, a low signal (asserted) on pin 5 causes the actuator to move to Position A and a high signal (deasserted) on pin 5 causes the actuator to move to Position B. However, in Mode 4 the actuator is normally enabled when pin 6 is in its high (deasserted) state. It is disabled only when pin 6 is pulled low (asserted), which is the opposite of how it works in Mode 3.

Mode Setup

To set the actuator mode, connect it to an RS-232 serial port as described in the section **Establishing Serial Communications**. To see the current setting, enter the **SM** command as shown in the **Serial Commands** chart on page 5. To change the mode, enter **SMn**, where n is a number between 1 and 4.

Mode settings are saved when the actuator is powered off.

SETTING THE MODEL SWITCH

The setting of the switch determines the serial port response seen to the **VR** command. Most end users will never need to perform this procedure, since the control module will be properly configured for the motor/gearbox assembly it shipped with. However, it is relevant for instrument builders who keep a stock of control modules which will be paired with a variety of motor/gearbox assemblies.

NOTE: If the switch is set to an invalid position (other than 1 through 5) the position indicator lights will flash continuously and the configuration will not change. If the unit is powered up with the model switch set in an invalid position, the VR command will elicit the response: **Not valid! Select 1-6!**

1. Find the tag on the motor/gearbox assembly which identifies it as a Model EQ, EH, EP, ED,ET, or EB.
2. Follow the instructions for removing the control module cover in "Setting the Serial Port Configuration Switch" on page 6.
3. Locate the 10-position rotary switch marked SW3 (**Figure 4**) and set it according to the legend printed underneath the switch for the motor/gearbox assembly to which it will be coupled.
4. Snap the cover back on the control module. You will probably wish to change the configuration tag on the bottom of the control module to reflect this modification.