




Universal Actuator BCD Connection to an Agilent 8890 GC

Cable Connections

1. Connect the VICI 8890 BCD cable (I-26476) to the Valco universal actuator and to the port labeled "BCD" on the 8890.
2. Plug in the power cord of the universal actuator.
5. Consult the Agilent 8890 user manual as required to configure valve #7 as a multiposition valve, with Invert BCD set, and a time delay of at least 2 seconds.

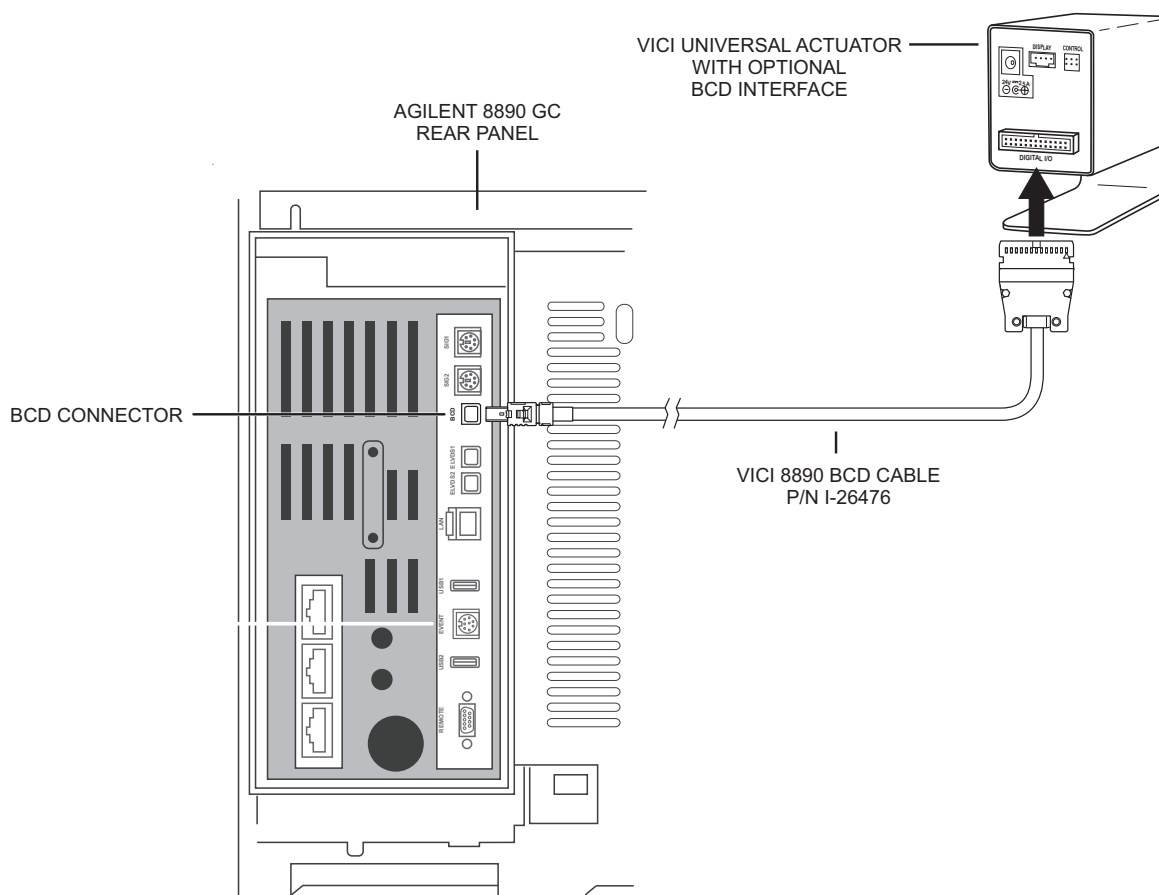

North America, South America, and Australia/Oceania contact:

VICI® Valco Instruments Co. Inc.
 P.O. Box 55603
 Houston, TX 77255
 Sales: (800) 367-8424
 Tech: (713) 688-9345
 Fax: (713) 688-8106 valco@vici.com

Europe, Asia, and Africa contact:

VICI® VICI AG International
 Parkstrasse 2
 CH-6214 Schenkon
 Switzerland
 Phone: +41 41 925 6200
 Fax: +41 41 925 6201 info@vici.ch

Cheminert® and VICI® are registered trademarks of Valco Instruments Co. Inc. and VICI AG